

Installing and Using our Cleanroom Tacky Mats




1 Ensure that the floor surface is clean, dust-free and dry before installation.




2 For a smoother and more secure mounting surface, we recommend using one of our [Stainless Steel Tacky Mat Frames](#).



3 Use the adhesive strips on the underside to secure frame to floor. You may also consider using mechanical fixings for additional security.

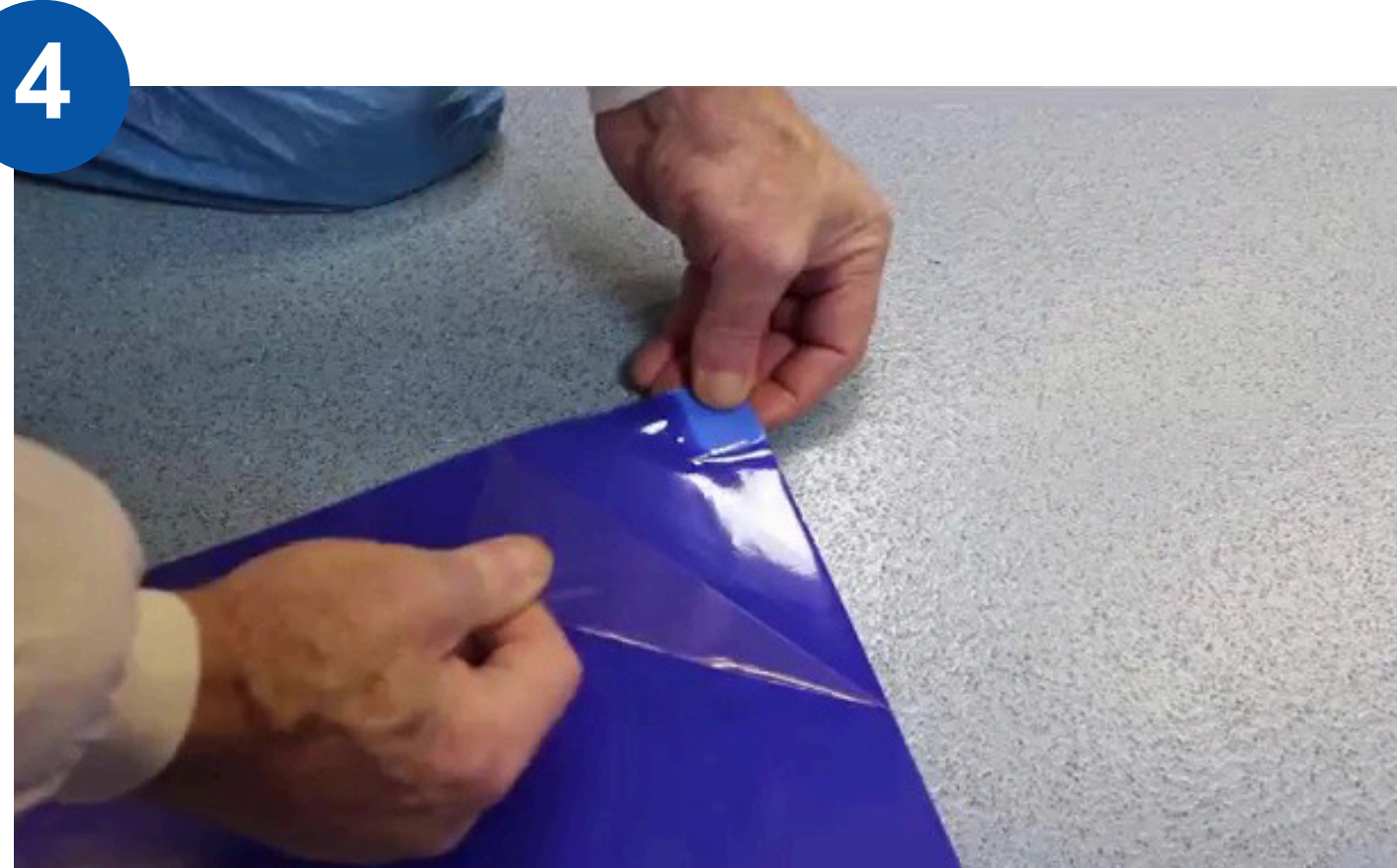


4 Peel back the protective film on the base of the tacky mat.

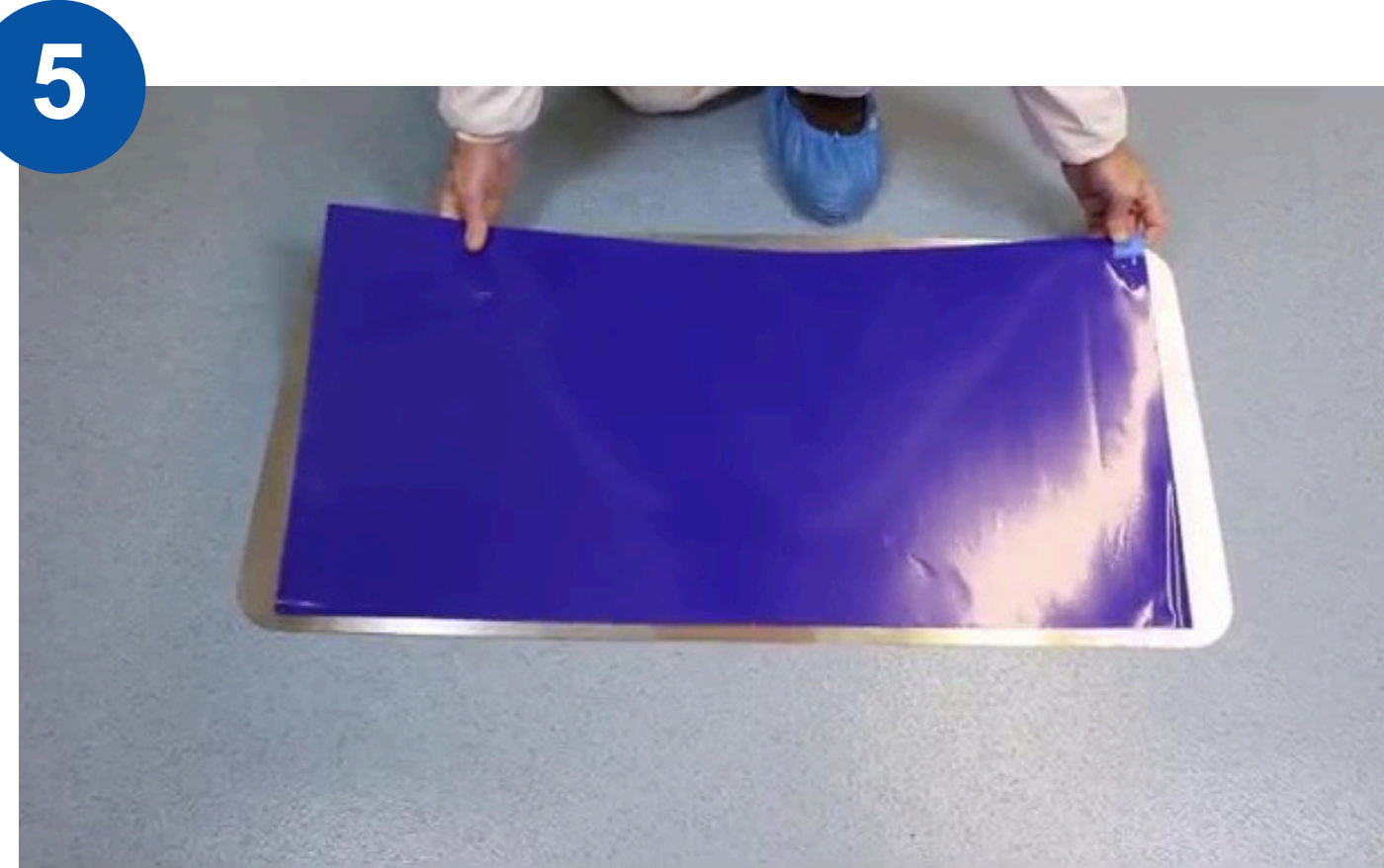


5 Place the mat on top of the frame in your desired location.




6 Before removing the top layer, use your hands to firmly smooth the tacky mat in place all over.

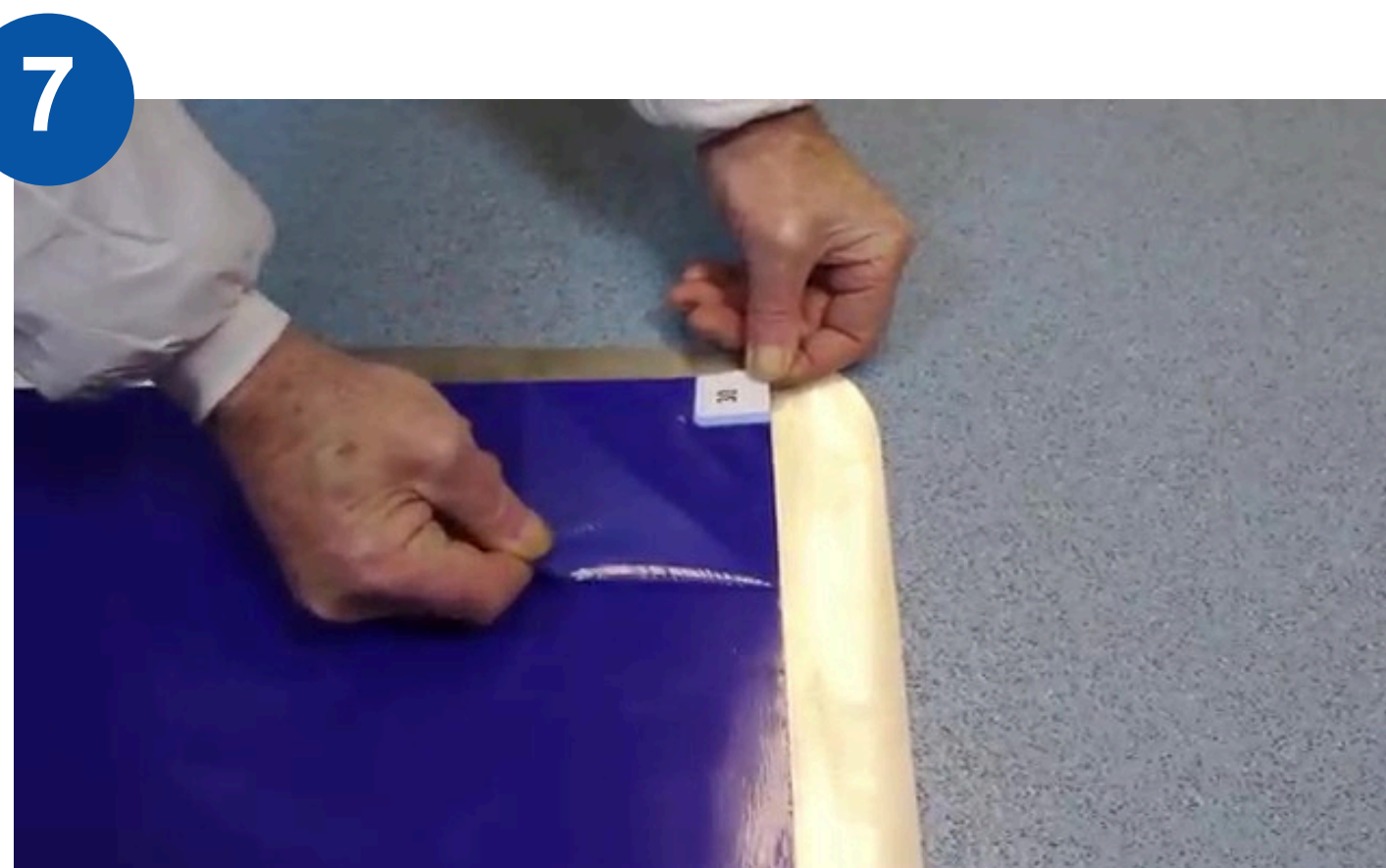


7 Peel back the protective top layer of the mat, being careful not to remove the first tacky layer.

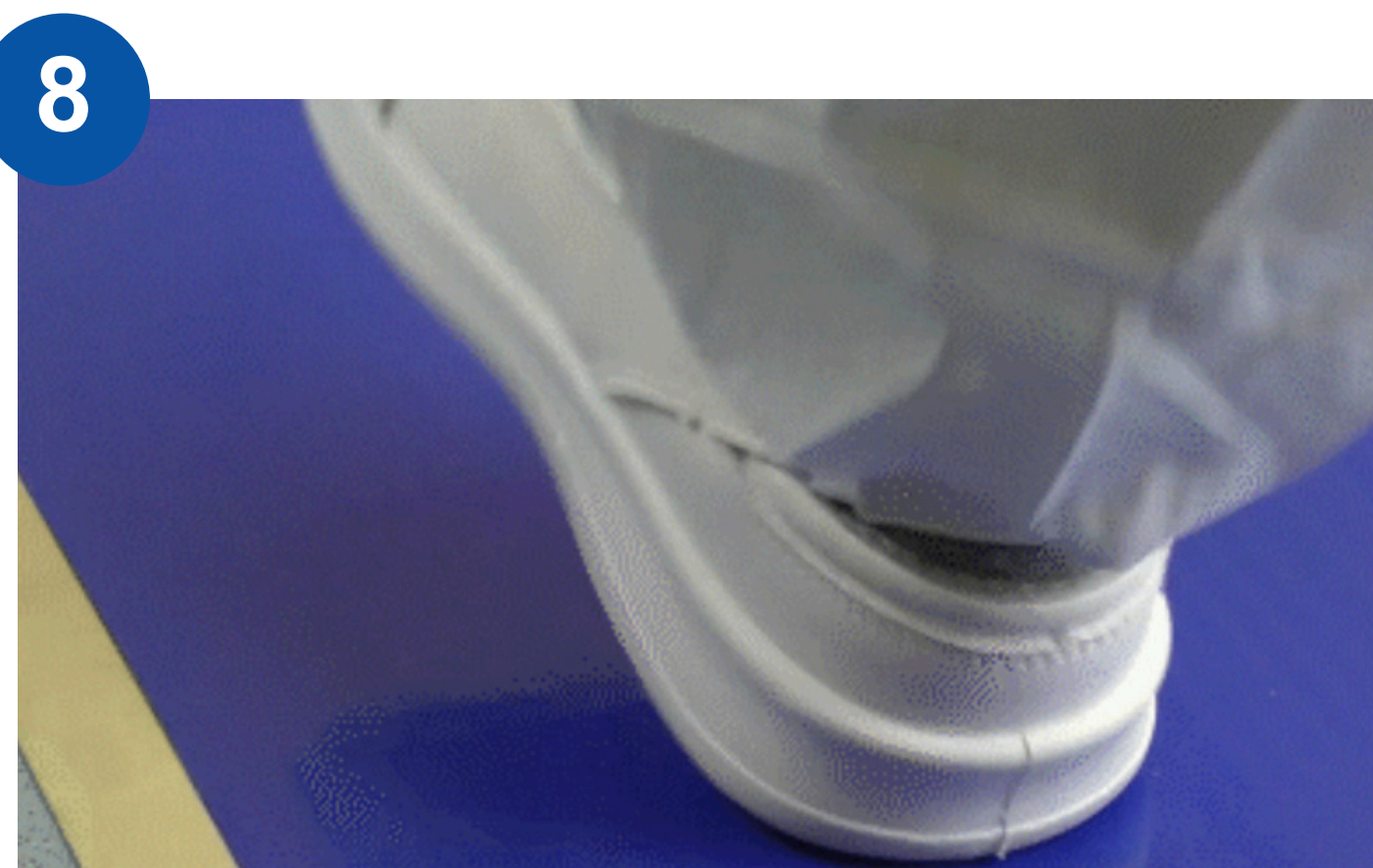

8 When walking across the mat, please take time to walk in a slower, deliberate manner.

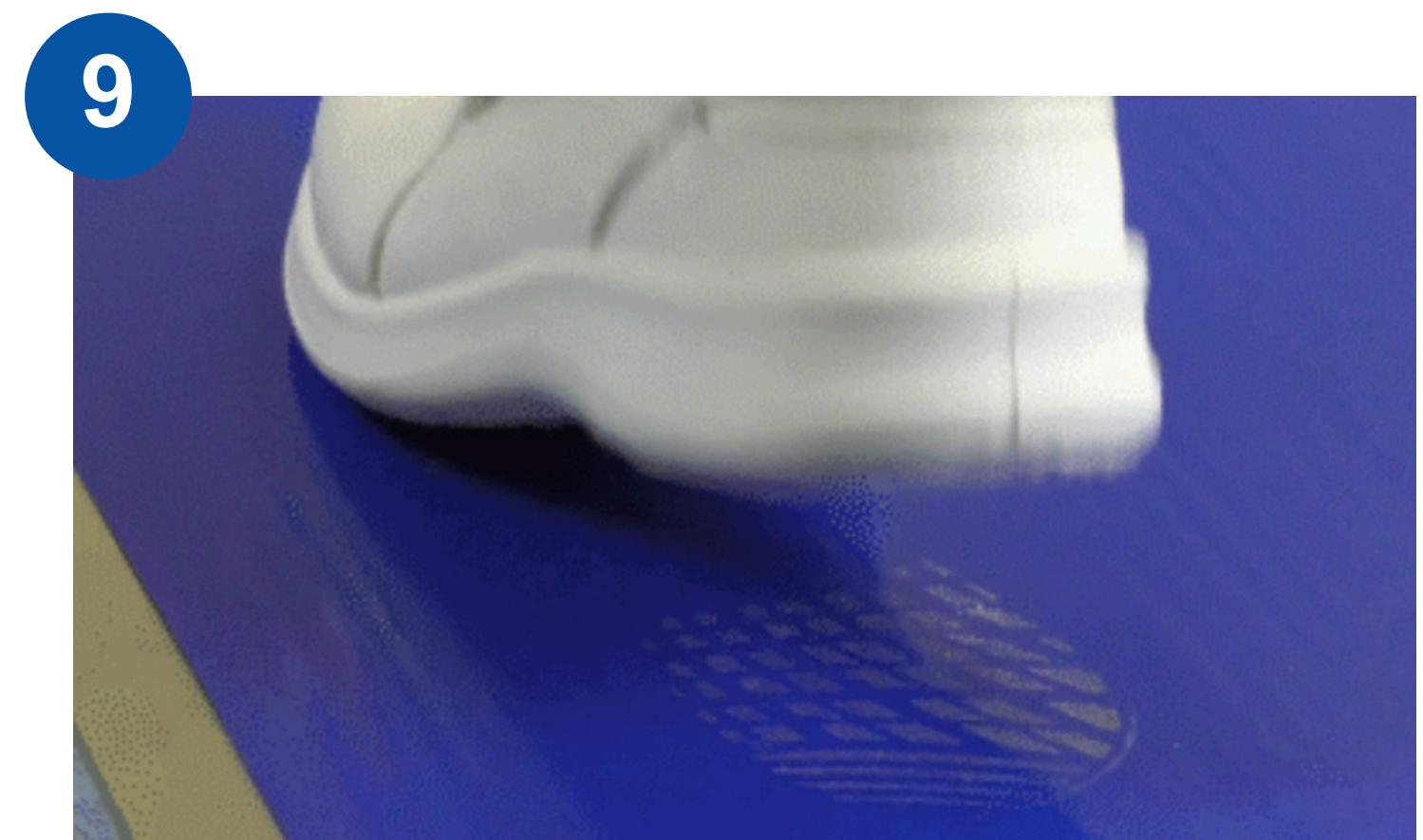

9 We recommend rocking your foot from heel to toe as you walk, pressing your foot firmly into the mat.

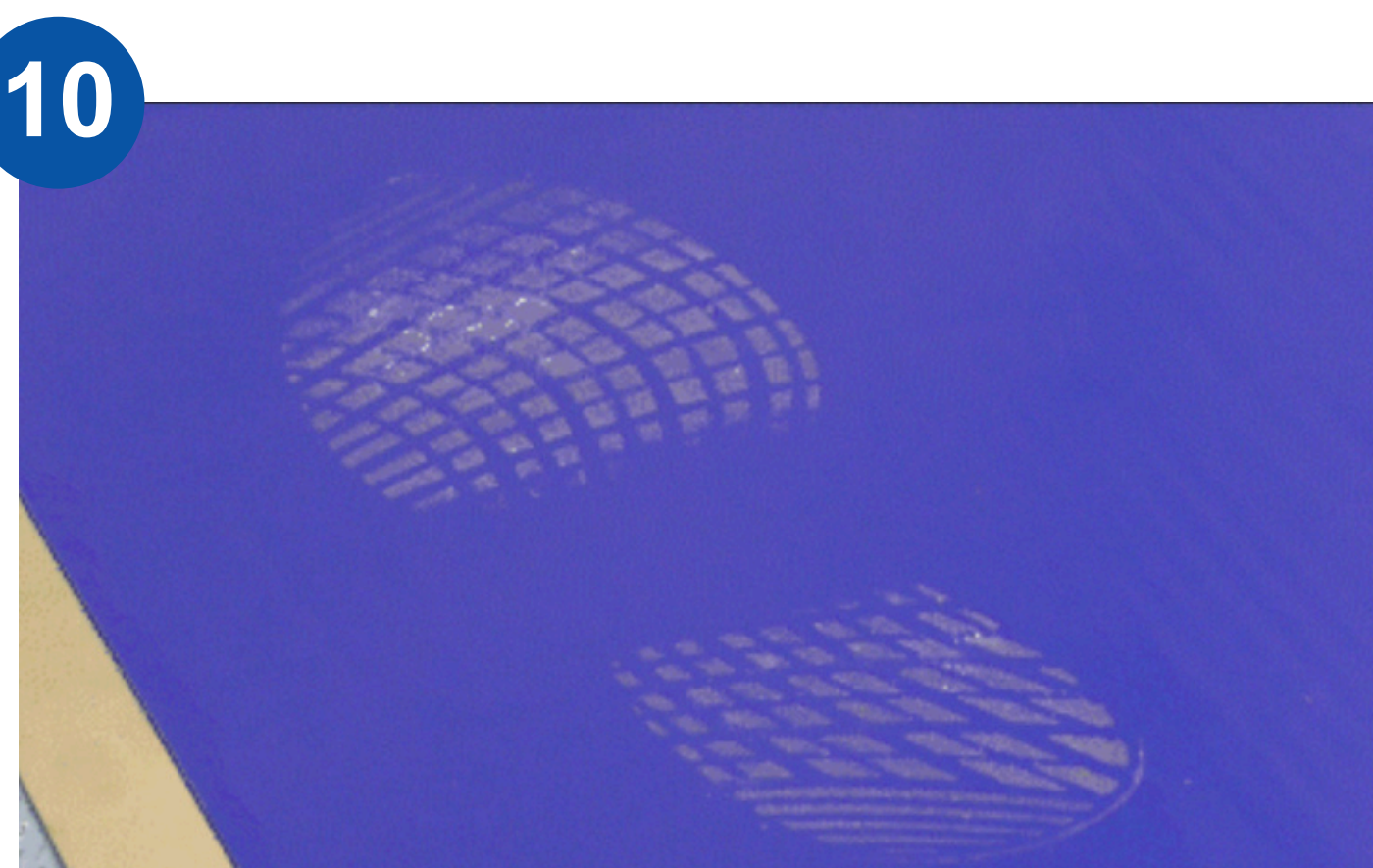

10 When the layer is visibly contaminated and losing its tack, you can remove the top layer from the numbered corner.