

Crystal

TITANIUM

The fastest acting sterile sporicidal disinfectant



User Guide

Crystel

TITANIUM

RECOMMENDED USE

Titanium is ideal for surface disinfection within controlled environments. It can be used on workbenches, shelves, equipment, printers, heat sealers and surfaces outside of the LAF cabinet for a daily deep clean and as part of a rotational regime.


CONTENTS

- 300ml Activator (1g/kg sodium chlorite)
- 300ml Base (50g/kg citric acid)

Titanium generates Chlorine dioxide (ClO_2).

Once empty, dispose of container and contents in accordance with local policy and national regulations.


HOW TO ASSEMBLE TITANIUM

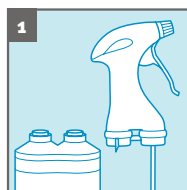


Take one Titanium pack.
Remove both bottle caps.



Place the head into the Titanium pack by inserting the tube into the open bottle neck.

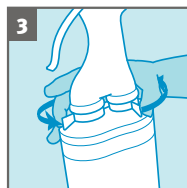

Attach the head to the Titanium pack by fastening the levers on each side.

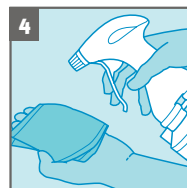

Titanium is now ready to use. It can either be applied directly onto a surface or onto a wipe.

MANUFACTURED BY

Tristel Solutions Limited. Cambs, CB8 7NY, UK

T +44 (0)1638 721500 **F** +44 (0) 1638 721911

E mail@tristel.com **W** www.tristel.com

Titanium is a good idea from **Tristel**